Operation Manual

Greiner Bio-One Mini Block Heater
REF 848050 (US) / 848060 (UK) / 848070 (EU)

Introduction:

Thank you for purchasing the Greiner Bio-One Mini Block Heater. Please read this manual thoroughly prior to operating the instrument. Using the device inconsistently with the instructions may result in damaging the device.

Text labeled with this sign contains safety notes. Read these safety precautions before using the device.

Intended Use:

The Mini Block Heater is intended only for indoor use and as a general laboratory tube heating device (dry bath). The digital block heater has an extremely compact design and provides excellent temperature uniformity and accuracy for heating of tubes from 0.2 mL up to 50 mL.

Specifications:

| | |
|-------------------|--------------------------|
| Temp. Range: | Ambient + 5 °C to 100 °C |
| Temp. Accuracy: | +/- 0.5 °C |
| Temp. Increments: | 0.1 °C |
| Temp. Uniformity: | +/- 0.2 °C |

| | |
|----------------------------------|----------------------------------------------------------------|
| Block Construction: | High Grade Aluminum |
| Operating Environment: | + 4 °C to + 45 °C + 39.2 °F to + 113 °F |
| Dimensions (Width/Depth/Height): | 4.4 (W) x 5.9 (D) x 4 (H) in. 11.2 (W) x 15 (D) x 10 (H) cm |
| Weight: | 2.0 lbs / 0.9 kg |
| Electrical: | 100 V to 240 V, 50-60 Hz |

Installation:

Place the Mini Block Heater on a flat and stable surface nearby an electrical outlet. It should not be near a heat or water source or in direct sunlight. Allow sufficient room around the unit for access and cooling. Six inches or 15 cm minimum on all sides is suggested. Before plugging in the Mini Block Heater compare your power supply with the electrical requirements.

Safety Precautions:

CAUTION: Hot surfaces, especially on the block, can cause serious injury or burns.

CAUTION: Do not put water or liquids into the well as shock, serious injury and death may occur.

CAUTION: Do not heat flammable or explosive substances as serious injury and death may occur.

Use of this product in any manner not specified by the manufacturer or modification of the product may cause injury and/or may void the warranty.

Do not use devices that have not been correctly installed or repaired (see also **Service and Contact**). Do not touch area around block or block well when unit is hot.

The Mini Block Heater should only be used for the specified applications. It must not be operated in hazardous, flammable or explosive environments and must not be used with explosive or highly reactive substances.

When handling toxic, radioactive or pathogenic liquids, observe the safety regulations of the country in question. If such liquids are spilled, the Mini Block Heater

must be cleaned carefully and properly after the unit has cooled down. Before using cleaning or decontamination methods others than those indicated in the manual, contact the manufacturer to ensure that the intended methods will not damage the device.

Tubes should be visually inspected for material damage prior to insertion. Damaged tubes must not be utilized. Make sure that lids of the tubes are tightly closed. When moving the Mini Block Heater from a cold room to a normal lab, allow it to warm up in the lab, but do not plug in the device in order to prevent damage caused by condensation.

Operation:

- Attach the power cord and place the power switch into the on “-” position.
- Remove the clear cover and install the appropriate block (purchased separately). Once installed, load your samples into the block and attach the cover.

Please note: Some blocks prevent the closing of the cover.

- Press the “SET” key to move the cursor and select the desired temperature using the up and down arrows.
- The instrument will now quickly heat to the desired temperature. The right side of the display will read “NOT OK” until the desired temperature has been reached. The display will then read “OK”.
- There is an over-temperature protection fuse that will cut heating in the event of an overheat. This means that once this fuse reaches temperatures above 125 °C, the heater will shut down until the temperature has decreased.

Cleaning and Maintenance:

Make sure that the dry bath and block are cool and the power cord is disconnected before performing any cleaning or maintenance. The outside of the Mini Block Heater can be cleaned with a moist cloth or 70 % isopropanol/water mixture. Do not immerse the Mini Block Heater in liquid or pour liquids over it. Excessive liquids should be avoided. The aluminum blocks can be removed for cleaning. Once removed, the blocks

can be cleaned with 70 % isopropanol/water mixture or can be sterilised in an autoclave at 250 °F / 121 °C for 20 minutes. After cleaning, ensure that all parts are dried thoroughly before attempting to reinsert in the unit.

Should a spill of toxic, radioactive or pathogenic liquids occur, the Mini Block Heater must be decontaminated thoroughly. Do not use any cleaning or decontamination methods except those recommended in this manual.

Accessory Blocks:

The following blocks are available for purchase; note that items marked with * are not compatible with the unit's cover:

| REF | Description |
|---------|--------------------------------------------------------------------------------------------------------------------------------|
| 848 902 | Block, 0.2 mL tubes or 5 PCR Strips of 8 tubes each |
| 848 913 | *Block, 4 mL / 5 mL Cryo.s™; 5 mL (12x75 mm), 4.5 mL (12x75 mm), 7 mL (13x100 mm) tubes |
| 848 915 | *Block, 15 mL conical tubes, 12 mL (16x100 mm), 12 mL (17x100 mm), 20 mL (16x152 mm), 14 mL (17x95 mm), 14 mL (18x95 mm) tubes |
| 848 916 | Block, 1.5 mL Reaction tubes, 0.5 mL PCR tubes |
| 848 921 | Block, 1 mL / 2 mL Cryo.s™; 4 mL (12x55 mm) tubes |
| 848 923 | Block, 1.5 mL / 2 mL Reaction tubes; 0.5 mL PCR tubes |
| 848 950 | *Block, 50 mL conical tubes |

Service and Contact:

This product is warranted to be free of defects in material and workmanship for a period of 2 years from the date of purchase. The warranty is valid only if the product is used in its intended purpose and within the guidelines specified in this instruction manual. In the event that service or technical support is required, please contact your nearest Greiner Bio-One office. When returning the Mini Block Heater, ensure that the device is fully decontaminated and does not present

any kind of health risk to our service personnel.

Disposal:

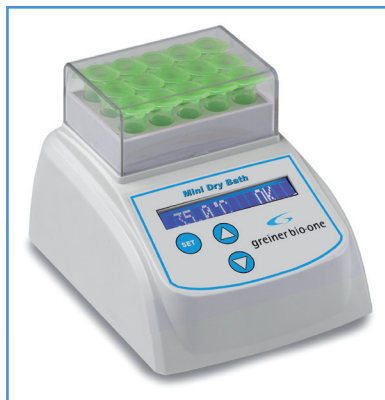
In the case the product is to be disposed of, the relevant legal regulations are to be observed. The disposal of electrical devices is regulated within the European Community by national regulations based on ED directive 2002/96/EC on waste electrical and electronic equipment (WEEE). As disposal regulations within EU may vary from country to country, please contact your supplier if necessary.

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Greiner Bio-One Co. Ltd.
Greiner Bio-One B.V.
Greiner Bio-One GmbH
Greiner Bio-One Ltd.
Greiner Bio-One North America Inc.