Operation Manual

Greiner Bio-One Mini Centrifuge
REF 843050 (US) / 843060 (UK) / 843070 (EU)

Introduction:

Thank you for purchasing the Greiner Bio-One Mini Centrifuge. Please read this manual thoroughly prior to operating the instrument. Using the device inconsistently with the instructions may result in damaging the device.

Text labeled with this sign contains safety notes. Read these safety precautions before using the device.

Intended Use:

The Mini Centrifuge is intended only for indoor use and for separating aqueous solutions and suspensions of various densities predominantl in research laboratories. It is designed for quick spin downs of approved microcentrifuge tubes (0.2 mL - 2.0 mL) as well as eight-position PCR strips.

Specifications:

Speed:	6000 rpm / 2000 x g
Capacity:	8 x 1.5/2.0 mL 32 x 0.2 mL or 4 x PCR strips
Dimensions (Width/Depth/Height):	4.5 (W) x 5.9 (D) x 4.5 (H) in. 11.4 (W) x 15 (D) x 11.4 (H) cm

Electrical:	100 V to 240 V, 50-60 Hz
Max. Radius:	1.93 in. / 4.9 cm
Weight:	2.2 lbs. / 1 kg
Operating Environ- ment:	+ 4 °C to 45 °C + 39.2 °F to 113 °F [max. 80 % relative humidity]

Installation:

Place the Mini Centrifuge on a stable, horizontal and level surface near an electrical outlet. In accordance with the safety regulations of IEC 1010-2-020 a safety distance of 30 cm should be observed around the centrifuge during operation. The Mini Centrifuge should not be near a heat or water source or in direct sunlight. Before plugging in the centrifuge, compare your power supply with the electrical requirements. Before starting the centrifuge for the first time, make sure the rotor is securely attached.

Safety Precautions:

The rotor must always be securely attached. If the centrifuge makes unusual noises when started, switch off the device immediately. The rotor must only be loaded symmetrically (see also section **Loading the Rotor**). Do not move the centrifuge during the run. Do not use centrifuges that have not been correctly installed or repaired (see also **Service and Contact**). The Mini Centrifuge should only be used for the specified applications. It must not be operated in hazardous or flammable environments and must not be used to centrifuge explosive or highly reactive substances.

When handling toxic, radioactive or pathogenic liquids, observe the safety regulations of the country in question. If such liquids are spilled, the centrifuge must be cleaned carefully and properly.

Before using cleaning or decontamination methods others than those indicated in the manual, contact the manufacturer to ensure that the intended methods will not damage the centrifuge.

Tubes should be visually inspected for material damage prior to centrifugation. Damaged tubes must

not be centrifuged. Make sure that lids of the tubes are tightly closed before centrifuging. When moving the centrifuge from a cold room to a normal lab, allow it to warm up in the lab, but do not plug in the centrifuge in order to prevent damage caused by condensation.

Rotor Exchange:

The Mini Centrifuge includes a second rotor that is designed for holding PCR tubes (0.2 mL) and PCR strips (8-position). The rotor is conveniently stored in a compartment on the bottom of the centrifuge (see picture). To access this rotor, turn the machine upside down and remove it from the compartment.

- **Removing the Rotor from the motor shaft:**
The new Clik-n-Lok™ rotor attachment system allows for instant exchange of rotors. To remove a rotor, simply pull it upwards, off the shaft. (NO TOOL IS REQUIRED.)
- **Attaching the Rotor to the motor shaft:**
Place the rotor on the motor shaft and press down until the rotor comes to a stop. The rotor is now secured. (NO TOOL IS REQUIRED)

Operation:

- Attach the power cord and place the power switch into the on “-” position.
- Open the lid and load your samples into the rotor. Always ensure a balanced load. Please see the following section, **Loading the Rotor**, for details.
- Close the lid. The rotor quickly accelerates to the maximum speed of 6000 rpm.
- When the desired run time is completed, push on the lid button to open the lid. The rotor quickly brakes to a stop and the samples can be retrieved.


Loading the Rotor:


To ensure safe operation and long life of the instrument, samples must always be loaded symmetrically into the rotor in a balanced fashion.

Tubes and strips must always be loaded symmetrically across from one another and contain the same volume of liquid. See below:

Example of properly balanced rotor: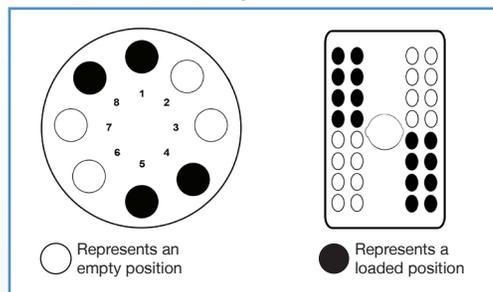

Example of improperly balanced rotor:

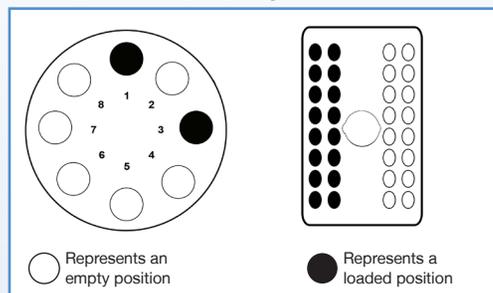Cleaning and Maintenance:

The outside of the Mini Centrifuge can be cleaned with a moist cloth and a mild noncorrosive detergent. Do not immerse the centrifuge in liquid or pour liquids over it. Excessive liquids should be avoided and liquids should not come in contact with the motor. Make sure that the centrifuge is disconnected from the main supply during cleaning. The centrifuge rotors can be removed for cleaning. Once removed, the rotors can be cleaned with 70 % isopropanol/water mixture or can be sterilised in an autoclave at 250 °F/121 °C for

20 minutes. After cleaning, ensure that all parts are dried thoroughly before attempting to operate the unit. Allow the rotor once to accelerate to the maximum speed. Should a spill of toxic, radioactive or pathogenic liquids occur within the rotor or the rotor chamber, the centrifuge must be decontaminated thoroughly. Do not use any cleaning or decontamination methods except those recommended in this manual.

Service and Contact:

This product is warranted to be free of defects in material and workmanship for a period of 2 years from the date of purchase. The warranty is valid only if the product is used in its intended purpose and within the guidelines specified in this instruction manual. In the event that service or technical support is required, please contact your nearest Greiner Bio-One office. When returning the centrifuge, ensure that the device is fully decontaminated and does not present any kind of health risk to our service personnel.

Tightening the Set Screw of the motor hub:

For self-tightening the set screw first remove the rotor from the motor shaft by simply pulling it upwards. If the rotor hub is loose, tighten the little screw on the side of the hub with a 1.5mm Allen wrench (make sure there is at least 2-3mm of space between the bottom of the hub and the base of the motor shaft). After snap on the rotor again. Make sure the unit is plugged in and turned On. Close the lid and ensure that the rotor begins to spin. Allow to spin for 10 seconds and ensure that the centrifuge does not show too much vibration and does not “walk” on the bench-top.

Disposal:

In the case the product is to be disposed of, the relevant legal regulations are to be observed. The disposal of electrical devices is regulated within the European Community by national regulations based on ED directive 2002/96/EC on waste electrical and electronic equipment (WEEE). As disposal regulations within EU may vary from country to country, please contact your supplier if necessary.

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