Operation Manual

Greiner Bio-One Mini Vortex Mixer
REF 845050 (US) / 845060 (UK) / 845070 (EU)

Introduction:

Thank you for purchasing the Greiner Bio-One Mini Vortex Mixer. Please read this manual thoroughly prior to operating the instrument. Using the device inconsistently with the instructions may result in damaging the device. The device should be maintained only at an authorised service center.

Text labeled with this sign contains safety notes. Read these safety precautions before using the device.

Intended Use:

The Mini Vortex Mixer is a compact, powerful, vortex mixer designed for vortexing tubes of various sizes (0.4 mL to 50 mL). The Mini Vortex Mixer is intended only for indoor use predominantly in research laboratories.

Specifications:

Speed:	Fixed: 2800 rpm
Mode:	TOUCH mix
Head:	Rubber cup standard for single tube mixing
Orbit:	4 mm / 0.157 in.
Dimensions (Width/Depth/Height):	9.4 x 9.9 x 6.6 cm 3.7 x 3.9 x 2.6 in.

Weight:	1 lbs. / 0.4 kg
Electrical:	100 to 240 V, 50-60 Hz
Operating environment:	+ 4 °C to + 45 °C + 39.2 °F to + 113 °F

Installation:

Place the Mini Vortex Mixer on a clean, stable and horizontal surface nearby an electrical outlet.

Press the Mini Vortex Mixer down slightly to allow the feet to adhere. The Mini Vortex Mixer should not be near a heat or water source or in direct sunlight. Before plugging in the Mini Vortex Mixer compare your power supply with the electrical requirements

Safety Precautions:

The Mini Vortex Mixer should only be used for the specified applications.

It must not be operated in hazardous, flammable or explosive environments and must not be used with explosive or highly reactive substances.

When handling toxic, radioactive or pathogenic liquids, observe the safety regulations of the country in question. If such liquids are spilled, the centrifuge must be cleaned carefully and properly.

Before using cleaning or decontamination methods others than those indicated in the manual, contact the manufacturer to ensure that the intended methods will not damage the device.

Tubes should be visually inspected for material damage prior to vortexing. Damaged tubes must not be utilized. Make sure that lids of the tubes are tightly closed. When moving the Mini Vortex Mixer from a cold room to a normal lab, allow it to warm up in the lab, but do not plug in the device in order to prevent damage caused by condensation. Do not use devices that have not been correctly installed or repaired (see also **Service and Contact**).

Operation:

Take the tube/vessel to be mixed, and pressed downward against the head to activate the vortex mixing ac-

tion. To get the desired mix, vary the angle of contact and pressure against the mixing head. When finished mixing, lift the tube upward.

Cleaning and Maintenance:

No routine maintenance is required other than to keep the unit clean. Cleaning can be done with a damp cloth. Avoid the use of solvents as they may attack the rubber cup head or the ABS plastic housing. Do not immerse the Mini Vortex Mixer in liquid and do not pour liquids over it. Excessive liquids should be avoided. After cleaning, ensure that all parts are dried completely. Do not subject the Mini Vortex Mixer to temperatures outside of the recommended operating range.

Should a spill of toxic, radioactive or pathogenic liquids occur, the Mini Vortex Mixer must be decontaminated thoroughly. Do not use any cleaning or decontamination methods except those recommended in this manual.

Service and Contact:

This product is warranted to be free of defects in material and workmanship for a period of 2 years from the date of purchase. The warranty is valid only if the product is used in its intended purpose and within the guidelines specified in this instruction manual. In the event that service or technical support is required, please contact your nearest Greiner Bio-One office. When returning the centrifuge, ensure that the device is fully decontaminated and does not present any kind of health risk to our service personnel.

Disposal:

In the case the product is to be disposed of, the relevant legal regulations are to be observed. The disposal of electrical devices is regulated within the European Community by national regulations based on ED directive 2002/96/EC on waste electrical and electronic equipment (WEEE). As disposal regulations within EU may vary from country to country, please contact your supplier if necessary.

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Greiner Bio-One Co. Ltd.
Greiner Bio-One B.V.
Greiner Bio-One GmbH
Greiner Bio-One Ltd.
Greiner Bio-One North America Inc.