# Operation Manual

Greiner Bio-One Vortex Mixer  
REF 844050 (US) / 844060 (UK) / 844070 (EU)

## Introduction:

Thank you for purchasing the Greiner Bio-One Vortex Mixer. Please read this manual thoroughly prior to operating the instrument. Using the device inconsistently with the instructions may result in damaging the device.

Text labeled with this sign contains safety notes. Read these safety precautions before using the device.

## Intended Use:

The Vortex Mixer is a compact, very powerful, vortex mixer designed for vortexing tubes of all various sizes (0.4 mL to 50 mL). It implements dynamic balance technology to provide smooth speed control and quiet operation with low excess vibration. The Vortex Mixer is intended only for indoor use predominantly in research laboratories.

## Specifications:

Speed:	User settable 200 to 3200 rpm
Mode:	ON (continuous), OFF and TOUCH mix
Head:	Rubber cup standard for single tube mixing

Feet:	Four no-walk suction cup
Orbit:	3 mm / 0.12 in.
Dimensions (Width/Depth/Height):	13 x 16 x 17 cm 5 x 6.3 x 6.7 in.
Weight:	8.4 lbs. / 3.8 kg
Electrical:	120 V/60 Hz, 230 V/50 Hz
Operating environment:	+ 4 °C to + 45 °C + 39.2 °F to + 113 °F

## Installation:

Place the Vortex Mixer on a clean, stable and horizontal surface nearby an electrical outlet.

Press the Vortex Mixer down slightly to allow the feet to adhere. The Vortex Mixer should not be near a heat or water source or in direct sunlight. Before plugging in the Vortex Mixer compare your power supply with the electrical requirements. Check that the cup head is securely fastened by gently rotating it clockwise.

Make sure that the mode switch is in the OFF position and plug the vortex mixer into a properly grounded outlet. Using the SPEED control knob, adjust the mixing speed to the desired setting. Start at the middle of the speed range if you are uncertain of which setting you need.

## Safety Precautions:

The Vortex Mixer should only be used for the specified applications.

It must not be operated in hazardous, flammable or explosive environments and must not be used with explosive or highly reactive substances.

When handling toxic, radioactive or pathogenic liquids, observe the safety regulations of the country in question. If such liquids are spilled, the centrifuge must be cleaned carefully and properly.

Before using cleaning or decontamination methods others than those indicated in the manual, contact the manufacturer to ensure that the intended methods will not damage the device.

Tubes should be visually inspected for material dama-

ge prior to vortexing. Damaged tubes must not be utilized. Make sure that lids of the tubes are tightly closed. When moving the Vortex Mixer from a cold room to a normal lab, allow it to warm up in the lab, but do not plug in the device in order to prevent damage caused by condensation. Do not use devices that have not been correctly installed or repaired (see also **Service and Contact**).

## Operation:

Set the mode switch to the desired setting of ON (continuous mixing) or TOUCH. In touch mode, the tube to be mixed is pressed downward against the head to activate the vortex mixing action. Touch the tube to be mixed to the head and observe the vortex mixing action. Adjust the SPEED knob to obtain more or less vortex mixing action.

## Head Removal and Replacement:

To change or replace the rubber head attachment, grasp the unit body with your left hand and then gently rotate the head attachment counter clockwise with your right hand until the cup head is completely unscrewed. Replace the head attachment by gently rotating clockwise until it is tight and you feel the rubber start to grip the top of the dynamic drive assembly.

## Cleaning and Maintenance:

No routine maintenance is required other than to keep the unit clean. Cleaning can be done with a damp cloth. Avoid the use of solvents as they may attack the rubber cup head or the ABS plastic housing. Do not immerse the Vortex Mixer in liquid and do not pour liquids over it. Excessive liquids should be avoided. After cleaning, ensure that all parts are dried completely. Do not subject the Vortex Mixer to temperatures outside of the recommended operating range.

Should a spill of toxic, radioactive or pathogenic liquids occur, the Vortex Mixer must be decontaminated thoroughly. Do not use any cleaning or decontamination methods except those recommended in this manual.

### Service and Contact:

This product is warranted to be free of defects in material and workmanship for a period of 2 years from the date of purchase. The warranty is valid only if the product is used in its intended purpose and within the guidelines specified in this instruction manual. In the event that service or technical support is required, please contact your nearest Greiner Bio-One office. When returning the centrifuge, ensure that the device is fully decontaminated and does not present any kind of health risk to our service personnel.

### Disposal:

In the case the product is to be disposed of, the relevant legal regulations are to be observed. The disposal of electrical devices is regulated within the European Community by national regulations based on ED directive 2002/96/EC on waste electrical and electronic equipment (WEEE). As disposal regulations within EU may vary from country to country, please contact your supplier if necessary.

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Greiner Bio-One Co. Ltd.
Greiner Bio-One B.V.
Greiner Bio-One GmbH
Greiner Bio-One Ltd.
Greiner Bio-One North America Inc.