

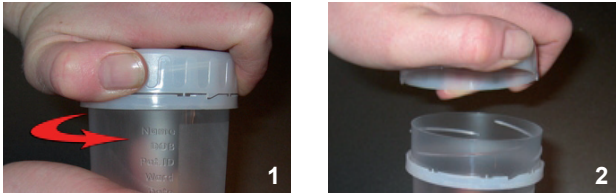

Polypropylene Multipurpose Beaker

with Integrity Seal

The Greiner Bio-One multipurpose beaker is manufactured from polypropylene. The beaker features an integrity seal, which ensures integrity and sterility of the interior. The beaker is leak-proof, unbreakable and resistant against many chemical influences.

The advantages of this hygienic all-rounder:

- Integrity seal guarantees sterility – it is no longer necessary to individually pack each beaker
- Time-saving, as there is no need to remove the beaker from a pack
- More information and better traceability – catalogue number, lot number and expiry date are printed on the base of the beaker
- With writing space and exact graduation marks

<p>Handling</p> <ol style="list-style-type: none"> 1. The lid is assembled on the beaker. To open, twist the lid thereby perforating the integrity seal. A light resistance must be overcome. 2. By opening the beaker, the sealing ring will be detached and the beaker is no longer sterile inside. 3. When closing the beaker with a new ring screw cap, a light resistance must be overcome. The pressure is necessary to ensure that the integrity seal of the screw cap snaps onto the beaker. This is a prerequisite for tight closure of the lid. If an audible “click” is heard, the closure has been applied correctly. 	
	

Ordering Information			
Cat. No.	Product Description	Quantity per Bag	Quantity per Case
724 461	Multipurpose Beaker with screw cap, 100 ml, graduated, natural, sterile	-	300
724 402	Multipurpose Beaker with screw cap, 100 ml, graduated, natural, non-sterile	-	300
724 401	Multipurpose Beaker bottom, 100 ml, without screw cap, graduated, non-sterile	-	720
724 170	Ring screw cap for Multipurpose Beaker cat. no. 724 401, natural, nonsterile	-	720