

CELLdisc™ Accessories

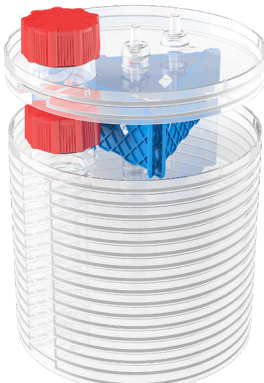

CELLlevator™ - Easy and secured CELLdisc™ stacking

The CELLlevator™ is a new CELLdisc™ accessory facilitating easy and secured stacking of two CELLdisc™ devices to enable best possible utilization of a given space (e.g. in an incubator).

The CELLlevator™ consists of three identical blue polypropylene segments which assembled form a uniform triangle. Centering ribs at the side of the segments recession guide the insertion assuring a simple and faultless assembly process. Thereafter the stacking device is positioned within the circumferential ring of the top plate of the CELLdisc™. Another CELLdisc™ can now be positioned on top of the CELLlevator™.

Ventilation of the lower CELLdisc™ is assured by the notches on the segments whereas planarity of the upper CELLdisc™ is assured by the elevations on the segments.

The maximum loading capacity of the assembled CELLlevator™ is 8 kg which corresponds a filled 40-layer CELLdisc™. Based on the application the CELLlevator™ is delivered unsterile but can be autoclaved if necessary.

<p>Key Facts</p> <ul style="list-style-type: none"> ➤ Easy and secured CELLdisc™ stacking ➤ Maximum loading capacity 8 kg ➤ Space saving storage ➤ Autoclavable (120 °C, 2 bar)* <p><small>*Autoclaving the CELLlevator™ more than 3 times is not recommended. After each autoclaving cycle the integrity of the segments has to be analyzed by the user.</small></p>		
--	---	--

Feature	Properties
Material	Polypropylene
Max. Loading Capacity	8 kg (refers to the assembled CELLlevator™)
Sterility	No
Autoclavability	Yes (120 °C, 2 bar), max. 3 times*
TP-range	4-37 °C

Ordering Information			
Order no.	Product Description	Qty. per bag	Qty. per case
878071	CELLlevator™ - Stacking device for CELLdisc™	1	9


For further information and/or sample ordering please visit our website www.gbo.com or contact us.

Germany (Main office): info@de.gbo.com | Austria: office@at.gbo.com | Belgium: info@be.gbo.com | Brazil: office@br.gbo.com | China: ts.bio@gbo-suns.com
 France: accueil@gbo.com | Hungary: office@hu.gbo.com | Italy: office@it.gbo.com | Japan: info@jp.gbo.com | Netherlands: info@nl.gbo.com
 Portugal: info@vacuette.pt | Spain: info@es.gbo.com | UK: info@uk.gbo.com | USA: info@us-gbo.com

CELLdisc™ Accessories

CELLring™ - Levelling ring compensating surface irregularities


CELLring™ assures an exact horizontal position of CELLdisc™ compensating surface irregularities, e.g. of a working bench or an incubator. Thus, cells and media distribute evenly in every single layer of your CELLdisc™.

<p>Key Facts</p> <ul style="list-style-type: none"> ➤ Ensures exact planar positioning of CELLdisc™ ➤ Guarantees consistent distribution of cells and media in every single layer 	
--	--

Ordering Information			
Order no.	Product Description	Qty. per bag	Qty. per case
878075	CELLring™ - Levelling ring compensating surface irregularities, e.g. of a working bench or an incubator	1	3

CELLhandle™ - Gripper for secure & easy transportation

CELLhandle™ facilitates secure and easy transportation of particularly large-sized CELLdisc™ formats as well as convenient emptying of these.

<p>Key Facts</p> <ul style="list-style-type: none"> ➤ Gripping device for easy lifting and emptying of large-sized CELLdisc™ ➤ Enables single-hand usage 	
---	--

Ordering Information			
Order no.	Product Description	Qty. per bag	Qty. per case
878074	CELLhandle™ - Gripper for secure and easy transportation	1	1

For further information and/or sample ordering please visit our website www.gbo.com or contact us.

Germany (Main office): info@de.gbo.com | Austria: office@at.gbo.com | Belgium: info@be.gbo.com | Brazil: office@br.gbo.com | China: ts.bio@gbo-suns.com
 France: accueil@gbo.com | Hungary: office@hu.gbo.com | Italy: office@it.gbo.com | Japan: info@jp.gbo.com | Netherlands: info@nl.gbo.com
 Portugal: info@vacuette.pt | Spain: info@es.gbo.com | UK: info@uk.gbo.com | USA: info@us-gbo.com